

24-Hour Emergency Number
(VelocityEHS) 1-800-255-3924 in U.S.

cleaning@alconox.com

alconox.com

- distributors
- technical information
- samples
- new developments
- white papers
- prices

DETOJET® Low-Foaming Liquid Detergent

- Concentrated to save you money
- Biodegradable and disposable
- Low-foaming heavy duty detergent for high performance results from your cleaning machinery
- Free rinsing to provide reliable results and no interfering residues
- For FDA cleaning validations and food contact surface cleaning

Used to clean: Hospital instruments, laboratory ware, optical components, pharmaceutical and cosmetics manufacturing equipment, tissue culture ware, plate glass, metal castings, forgings and stampings, industrial parts, animal cages, electronic components, tanks and reactors.

Used to remove: Soils, grit, grime, slime, grease, tars, resins, fats, oils, blood, tissue, particulates, inorganic residues, chemicals and solvents.

Surfaces cleaned: Recommended for glass, stainless steel, ceramic, cement and alkali-resistant plastics and metals such as stainless steel. Corrosion testing is advisable.

Cleaning method: Labware washers, dishwashers, washer-sterilizers, washer-sanitizers, cart washers, spray cabinet washers, electrocleaning, ultrasonic cleaners clean-in-place and manual cleaning.

Directions: Follow machine manufacturers directions. Not intended for powder detergent receptacles. If available, use liquid detergent reservoir, connecting tubes or metering pump. In general, use 1 – 2% solution in hot water wash cycle (3 – 5 Tbsp/gal, 1 1/4 – 2 1/2 oz/gal or 10 – 20 ml/L). Most machines need 2 1/2 oz added at a typical 2 gal. wash cycle. For soak, circulate or ultrasonic use, make a 1% solution in cold, warm or hot water. For difficult soils, raise water temperature and use more detergent. Wear protective gloves and eyewear when handling.

RINSE THOROUGHLY—preferably with running water. For critical cleaning, do final or all rinsing in distilled, deionized or purified water. For food contact surfaces, rinse with potable water. Used on a range of alkali resistant glass, ceramic, plastic and metal surfaces. Corrosion testing is advisable. Material compatibility guidelines available on request.

Convenient Sizes	Alconox Cat. #
Case of 4 x 1 gal (4 x 3.8L)	1601
5 gal (19L) jerrycan	1605
15 gal (57L) drum	1615
55 gal (208L) drum	1655
1 gal makes 200 gal (1 L makes 200L) solution	



DETOJET passes inhibitory residue test for water analysis and is FDA certified. DETOJET is available from leading laboratory, hospital, clinical and industrial suppliers. To find a distributor for Alconox Inc. detergents, use the Locate Dealers search tool at alconox.com. To request samples, use the Order Samples form at alconox.com. Or call Alconox Inc. at +914.948.4040, today.

DJTB.5

24-Hour Emergency Number
(VelocityEHS) 1-800-255-3924 in U.S.

cleaning@alconox.com

alconox.com

- distributors
- technical information
- samples
- new developments
- white papers
- prices

Physical Data	Typical Value
pH (as is)	12.4
Surface Tension 1% Sol'n (dyne/cm)	55
Conductivity 1% Sol'n (mS/cm at 21 degrees C)	0.7
Form	Liquid
Color	Clear
Solubility in Water	Completely soluble in all proportions
Hard Water Effectiveness	Highly effective
Biodegradability	Yes, meets EC648/2004
Foam Tendency	Low foaming
Shelf Life	42 months from end of month of manufacture
Fragrance and Dye Content	0%
Phosphate Content (as phosphorus)	10.1%
Flash Point (degrees F)	None
Density (lb/gal)	10.8
Organic Carbon (% calculated w/w)	0.5%

Chemical Description

Detojet is a highly alkaline detergent containing potassium hydroxide and silicate of soda. It should be used with caution and the precautions on the label should be followed. Detojet is surfactant free. Due to other components in the detergent mixture, only trace amounts of hypochlorite are present post production.

Cleaning Validation and Residue Detection

To quantify trace levels of Detojet for cleaning validation purposes, use conductivity, phosphate analysis, HPLC, IC, or TOC. For details see the Cleaning Validation page in Support at alconox.com. For lab accreditation tests to show the detergent is rinsed away, see the Accreditation Residue Detection in the Support section of alconox.com.

Health and Safety Information

OSHA Hazardous Ingredients	None
RCRA Hazard Class	Nonhazardous. Undiluted material may be hazardous and require regulatory compliance.
EPA Priority Pollutants	None
DOT Hazard Class	UN 1760, Corrosive Liquid, n.o.s., 8, PG II
Flammability	Nonflammable
Latex Content	None in product or packaging.
Irritation	Cat. 2 skin, Cat. 2A eye irritant
Oral Toxicity	LD ₅₀ > 500 mg/kg oral rat.
Inhalation Toxicity	Nonirritating
VOC Content	0%
GMO Content	None
BSE/TSE Source	None
2007/68/EC Annex IIIa Allergens	None
REACH Art. 57 (SVHC)	None
RoHS	No RoHS hazardous ingredients.
Carcinogenicity	NTP = No IARC = No OSHA = No
Ingredients on TSCA Inventory	Yes
EC 907/2006	Contains no enzymes, optical brighteners, perfumes, allergenic fragrances, or preservative agents.

Precautions

Avoid contact with skin, eyes and clothing. Do not take internally (see Directions). In case of contact immediately flush skin with plenty of water. For eyes, flush with plenty of water for at least 15 minutes and get medical attention immediately. A Safety Data Sheet is available at alconox.com.

Contact Alconox Inc. for purchase specifications. Information presented is typical. Not to be taken as specifications. Typical data is not a specification.

While the information in this report should not be considered to be a product warranty, we urge you to investigate, test and verify the suitability of Alconox Inc. detergents for your specific application. We, of course, can not give permission to use, or recommend the use of, our detergents where they infringe patents. No representation or warranty is made as to the safety of products or materials mentioned under the Federal Food Additives Amendment of 1958.

DJTB.5